DURAN®

Laborglaskatalog / Laboratory glassware Catalogue /
Catalogue de verrerie de laboratoire /
Catalogo de material para laboratorio

SCHOTT
glass made of ideas

Laborglas Laboratory Glassware Verrerie à emploi général Artículos de vidrio para laboratorio	1 7-22
Flaschen Bottles Flacons Frascos	2 23-36
Exsikkatoren Desiccators Desiccateurs Desecadores	3 37-44
Gläser für die Mikrobiologie Glassware for Microbiology Verrerie pour la microbiologie Vidrio para microbiología	4 45-56
Volumenmeßgeräte Volumetric Glassware Appareillage en verre pour la volumétrie Material volumétrico	5 57-68
Glasfiltergeräte, Trichter Glass Filtration Apparatus, Funnels Appareillage de filtration en verre, Entonnoirs Aparatos de vidrio de filtración, Embudos	6 69-90
Schliffbauteile Interchangeable Glassware Éléments interchangeables en verre rodé Piezas esmeriladas intercambiables	7 91-120
Weiterverarbeiterartikel Products for manipulators Articles et ébauches pour souffleurs de verre Artículos de procesamiento adicional	8 121-164

Becher, niedrige Form

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Beaker, low form

with graduation and spout
DIN 12331, ISO 3819

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Vasos, forma baja

graduados, con pico
DIN 12331, ISO 3819

Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	Originalpckg./Stück Stand. pack/quantity
21 106 07 ¹	5	22	30	10
21 106 08 ¹	10	26	35	10
21 106 14	25	34	50	10
21 106 17	50	42	60	10
21 106 24	100	50	70	10
21 106 29	150	60	80	10
21 106 36	250	70	95	10
21 106 41	400	80	110	10
21 106 48	600	90	125	10
21 106 53	800	100	135	10
21 106 54	1 000	105	145	10
21 106 63	2 000	132	185	10
21 106 68	3 000	152	210	4
21 106 73	5 000	170	270	1
21 106 86 ^{1, 2}	10 000	217	350	1

¹ Ohne Teilung

¹ Without graduation

¹ Sans graduation

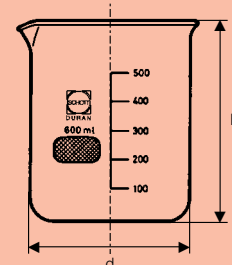
¹ Sin graduar

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² Non-DIN, non-ISO size

² Non conforme aux dimensions DIN et ISO

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Becher, hohe Form

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Beaker, tall form

with graduation and spout
DIN 12331, ISO 3819

Béchers, forme haute

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Vasos, forma alta

graduados, con pico
DIN 12331, ISO 3819

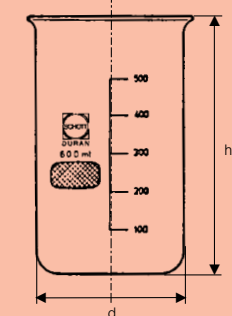
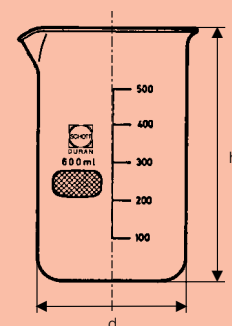
Best.-Nr. Cat. No.	Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	Originalpckg./Stück Stand. pack/quantity
21 116 17	21 117 17 ¹	50	38	70	10
21 116 24	21 117 24 ¹	100	48	80	10
21 116 29	21 117 29 ¹	150	54	95	10
21 116 36	21 117 36 ¹	250	60	120	10
21 116 41	21 117 41 ¹	400	70	130	10
21 116 48	21 117 48 ¹	600	80	150	10
21 116 53		800	90	175	10
21 116 54	21 117 54 ¹	1 000	95	180	10
21 116 63		2 000	120	240	10
21 116 68		3 000	135	280	2

¹ ohne Ausguß

¹ without spout

¹ sans bec

¹ sin pico

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CERAN® Protective laboratory plates

Saving in energy and time

The good permeability of CERAN® protective plates for infrared radiation results in heat energy transfer to the material to be heated with low losses.

This saves 20% and more time and energy. In addition, there is room for several items on the flat square area.

Chemically resistant

When working in the laboratory, boil-ing over and spilling of aggressive media is unavoidable in practice.

Even highly aggressive media can do no damage to the CERAN® lab-oratory protective plate.

Problem-free cleaning

The smooth pore-free surface of the CERAN® plate can be cleaned by hand and by machine without prob-

lems, and even in the same machine where your laboratory glassware is cleaned: in the dishwashing machine.

High temperature resistance

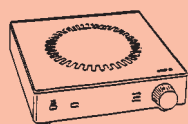
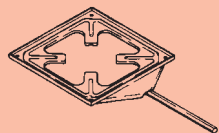
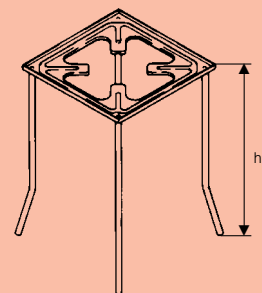
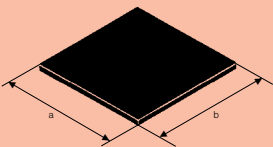
Application range from
– 200 °C to + 700 °C.
A particular advantage of CERAN® laboratory protective plates is their high continuous operation temperature.

Durability at
700 °C – 6,000 h
750 °C – 750 h
800 °C – 100 h

Even when the hot plate is quenched with cold water, there is no risk of fracture, since the resistance to temperature shock is greater than 650 °C. In order to avoid overheating, care must be taken that the above-mentioned limits are not exceeded when working with a bunsen burner. Our CERAN® protective laboratory plate retains its shape and flatness and dos not age.

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Quadrupods for CERAN® laboratory protection plates of chrome-nickel steel

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Plate holder for CERAN® laboratory protection plates of chrome-nickel steel

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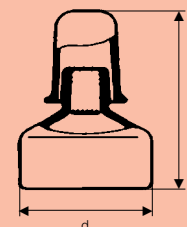
Spirit lamps, made of soda-lime

without filler tubulature
with ground-on cap

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Spirit lamps, made of soda-lime

with filler tubulature and stopper
NS 10, with ground-on cap

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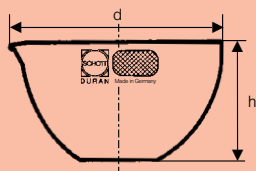
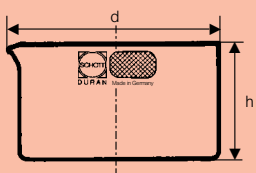
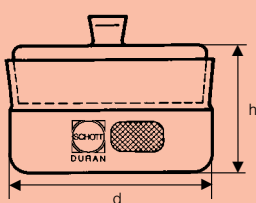
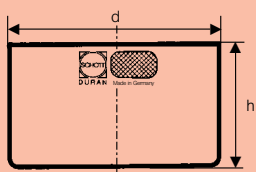
Cristallisoirs

Cristalizadores

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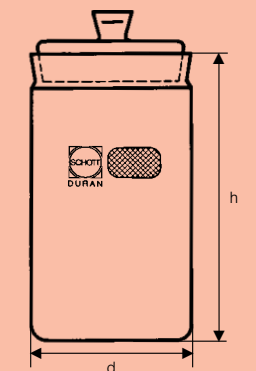
Weighing bottles with grip stopper, DIN 12605

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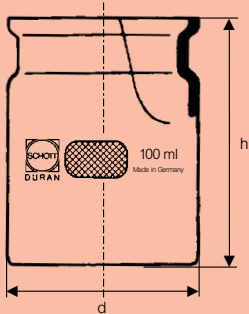
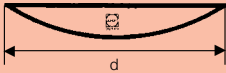
Watch glass dishes
with fused edges, DIN 12341

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Organ storage jars
without stoppers

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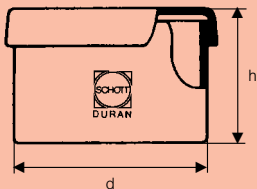
Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	Originalpkg./Stück Stand. pack/quantity
24 204 23	75	50	70	10
24 204 24	100	54	75	10
24 204 36¹	250	74	102	10

¹ Nur solange Vorrat reicht
¹ Discontinued, available until stocks exhausted
¹ Uniquement jusqu'à épuisement du stock
¹ Sólo hasta agotar existencias

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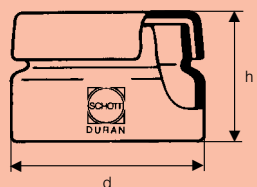
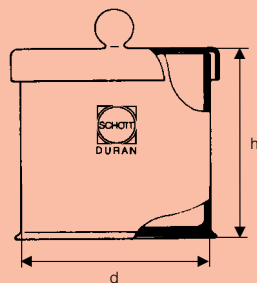
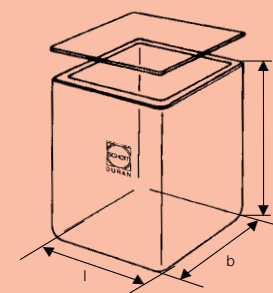
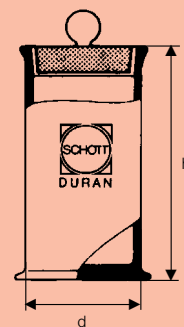
Cylinders with knobbed lid, polished rim

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¹ Nur solange Vorrat reicht

¹ Discontinued, available until stocks exhausted

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Spare cover plates upon request (minimum order: 10 pcs.).

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DURAN test tubes have a heavier wall. They therefore have considerable mechanical strength and excellent thermal shock resistance.

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SUPREMAX test tubes withstand temperatures up to 700 °C over prolonged periods, and even higher temperatures over short periods.

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FIOLAX test tubes have a thinner wall. They are therefore relatively resistant to rapid temperature change and local heating.

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mit Teilung

Laboratory bottles with DIN thread, ISO 4796
graduated

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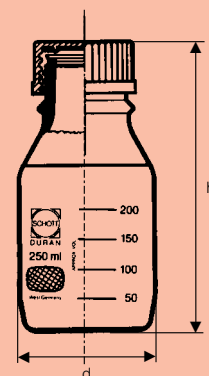
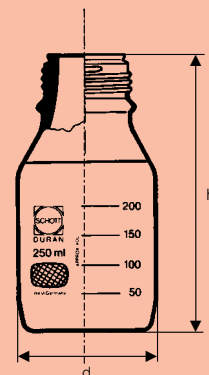
mit Teilung

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Laboratory bottles with DIN thread, ISO 4796, amber
graduated
without cap and pouring ring

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¹ A bead is moulded in the glass on the shoulder of the 100-2000 ml sizes to indicate the nominal capacity level.

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² bottle with **specially formed glass lip** for better pouring (this makes an additional plastic pouring ring **unnecessary**).

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Chemical and thermal properties see page 180.

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DURAN® laboratory bottles

DURAN® laboratory bottles are chemically resistant and stable. Supplemented by a plastic pouring ring they allow drip-free operations. Since all bottles from 100 ml onwards have only one size of screw thread, screw caps and pouring rings can be exchanged for one another. Bottle, pouring ring and screw top can be sterilised.

Properties:

Light protection

- by brown bottles up to 500 nm
- by plastic coating up to 380 nm
- usable as safe for chemicals

kept sterile by a safe seal for example

- of dairy products
- of serums

high resistance to temperature shock

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Further processing

The DURAN® laboratory bottles are manufactured from the tested borosilicate glass 3.3 and are suitable, for example, for fitting

- Screw thread pipes
- Connectors
- Tubes
- Sidearms.

Handling hints:

Freezing of substances

Recommendation: freeze the bottle in an inclined position (about 45 °) and only fill to at most 3/4 (to increase the surface area).

Temperature limit: – 40 °C, since the plastic caps and pouring rings do not withstand lower temperatures.

Melting of frozen substances

The frozen material can be melted by immersing the bottle into a liquid bath (temperature difference not more than 100 °C). In this manner, the frozen material is heated uniformly from every side without damaging the bottle. However, it can also be

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melted slowly from the top, so that the surface becomes liquid first and the material can expand.

Sterilisation

Important: during sterilisation, the screw cap may only be loosely (not more than one turn) screwed up (do not close fully). In a sealed bottle, no pressure release can take place. The resulting difference in pressure can result in fracture of the glass article (it will burst).

Further important notes:

Load dishwashing machines in such a manner that the glass articles – in particular the threads – do not bump into each other.

Pressure resistance:

DURAN® laboratory bottles are not suitable for vacuum and pressure works, except pressure resistant bottle **21 810 54**.

Information about pressure-resistant laboratory bottles can be supplied on request.

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Laboratory bottles, with DIN thread, ISO 4796, plastic coated (DURAN® Protect)²

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¹ A bead is moulded in the glass on the shoulder of the 100–2000 ml size to indicate the nominal capacity level.

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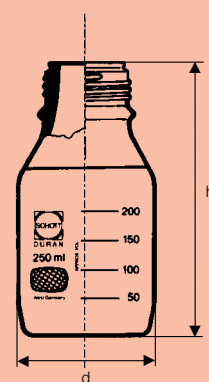
For spare-caps and pouring ring see page 27.

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Laboratory bottles with DIN thread

ISO 4796, **pressure resistant – 1/+ 1,5 bar**, blue graduation without cap and pouring ring

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¹ A bead is moulded in the glass on the shoulder to indicate the nominal capacity level.

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DURAN® Protect DURAN® laboratory bottles with plastic coating

Important notes:

The plastic coating provides protection against mechanical damage and it prevents leakage of the contents should the glass break.

The copolymer based plastic material remains transparent and adheres to the glass surface even after repeated sterilisation in an autoclave. (Steam

In the case of plastic-jacketed laboratory bottles, the autoclave instruction has to be observed when autoclaving with steam.

Autoclave instruction:

1. Autoclaving in steam max. 135 °C.
2. Switch off heating and let cool down slowly (20 min).
3. Aerate autoclave at 100 °C and remove water. Do not cool with distilled water or vacuum.
4. Open autoclave at 80 °C.

Microwave: plastic-coated laboratory bottles are not suitable for operations in a microwave oven.

Sterilisation

Important: during sterilisation, the screw cap may only be loosely (not

up to 135 °C; do not use distilled water for cooling.)

The special composition of the plastic material is largely UV absorbent (to 380 nm, s. page 173 transmission curves, Optical properties).

Pressure resistance

The bottles are not suitable for operations involving pressure or vacuum, except art. no. **21 815 54**.

Important:

Plastic-coated bottles are not suitable for freezing.

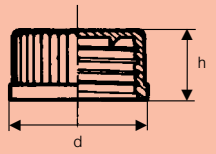
more than one turn) screwed up (do not close fully). In a sealed bottle, no pressure release can take place. The resulting difference in pressure can result in fracture of the glass article (it will burst).

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PP¹ Screw-caps, blue

with lip seal autoclavable up to 140 °C steam

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¹ Chemical and thermal properties see page 180.

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Bottles, reagent, wide neck
made of **soda-lime glass**,
DIN 12 038

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Bottles, reagent, wide neck
made of **soda-lime glass**,
DIN 12 039,
with standard ground joint

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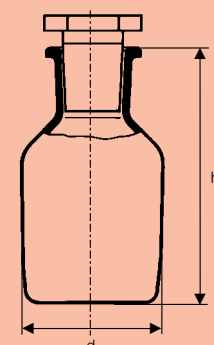
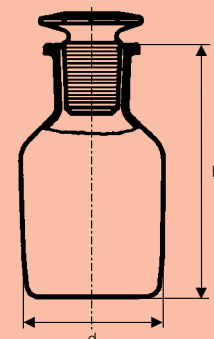
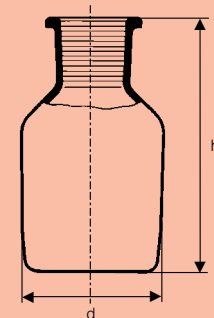
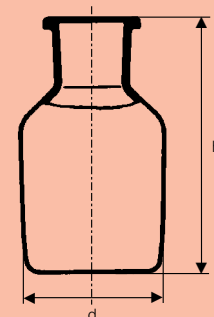
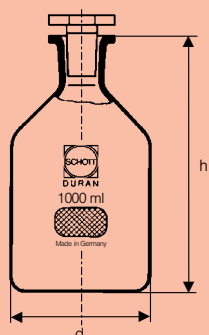
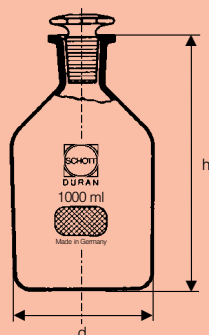
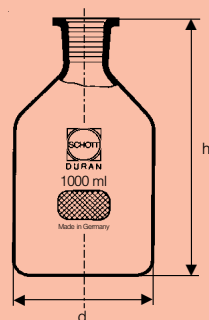
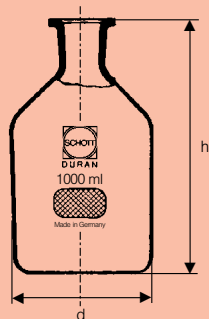
Important note: When DURAN glass stoppers are used with **soda-lime glass** reagent bottles, they may become stuck in the neck if temperature differences exceed 30 °C.

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² Bottles reagent, narrow neck, with neck 29/32, 1000 ml, can be used as burette bottle (reservoir bottle) see page 63.

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**DURAN bottles, reagent,
narrow neck**
DIN 12 036, ISO 4796
with standard ground joint

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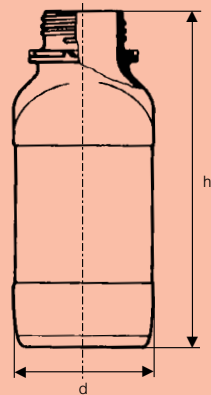
Square bottles,
narrow neck, with DIN thread, made
of **soda-lime glass, complete**
with dust cap and pouring ring

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Square bottles,
wide neck, with DIN thread,
made of **soda-lime glass,**
without cap and pouring ring

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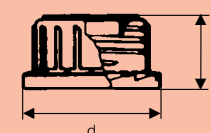
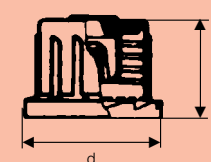
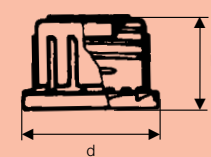
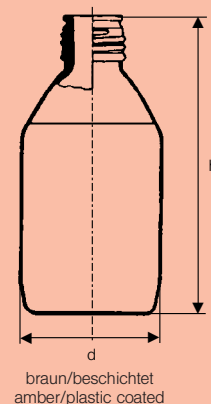
Round bottle, narrow neck,

amber/with and without plastic coated, with DIN-thread, made of **soda-lime glass, complete** with cap and pouring ring

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Tamper-proof PP screw-caps,

blue, for **narrow neck** square bottles

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Safety, tamper-proof PP

screw-caps, red for **narrow neck** square bottles

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Acid bottles, complete

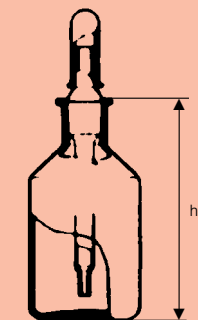
with NS-pennyhead stopper,
conicalshouldered,
interchangeable cap

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Dropping bottles

with interchangeable pipette NS,
soda-lime glass, complete with
rubber teats

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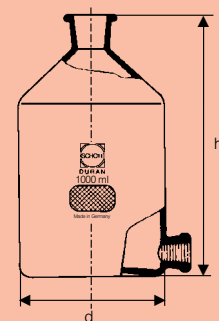
DURAN-Aspirator bottles,

NS tubulated, DIN 12037,
complete with NS stopcock
and NS stoppers

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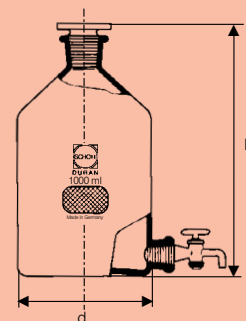
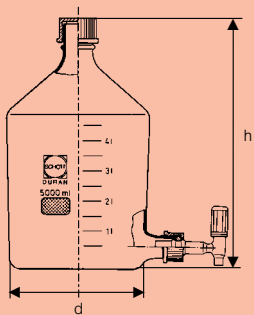
DURAN Aspirator bottles

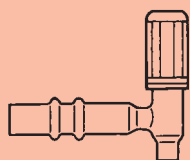
neck with GL 45 DIN screwthread, complete with GL 32 tubulature, screwthread adapter coupling, silicone gasket and draincock with PTFE spindle

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Draincock

with PTFE spindle for GL 32 DIN screwthread for aspirator bottle

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Desiccators

flange with groove and NOVUS
NS junction. Tubulature in lid,
with O-ring (PBR) and stopcock

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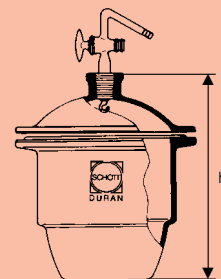
Desiccators

with flat flange and NOVUS NS
junction. Outlet in lid, with stopcock,
DIN 12 491

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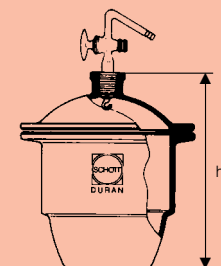
Desiccators

Delivery: base with NS side outlet
with stopcock and lid

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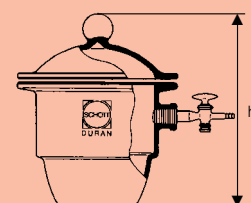
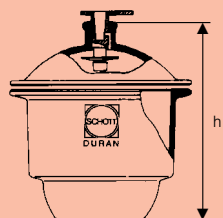
Desiccators

with **WERTEX** safety junction in lid,
DIN 12 491

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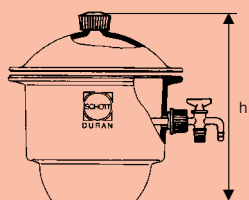
Desiccator accessories, such as closures and seals, see page 42/43.

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Desiccators

with **MOBILEX** screw thread
in lid and base, with stopcock

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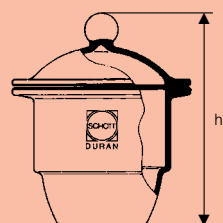
Desiccators

with plane flange, without junction,
with knobbed lid

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DURAN® desiccators – Variations and safety with DURAN

The range consists of five different desiccator types as well as all the appropriate spare parts such as bases, lids and a variety of accessories.

All desiccators and individual parts are made from the tried and tested DURAN borosilicate glass 3.3. Highest possible chemical resistance, minimum heat expansion as well as the consequent high resistance to temperature changes are among the characteristics of DURAN. These physical and chemical properties make DURAN ideal for use in laboratory glassware. You will find more information on DURAN borosilicate glass from SCHOTT on page 166 onwards.

DURAN desiccators come in two different versions – clear and plastic coated. The plastic coating of the DURAN Protect series is an additional protection for DURAN desiccators. It protects from scratches, mechanical damage and increases the stability of the item.

Five desiccator types are available, according to actual requirements:

- Desiccators with flat flange and no joint
- Desiccators with ring groove flange and NOVUS ground joint tubulature
- Desiccators with flat flange and NOVUS ground joint tubulature
- Desiccators with WERTEX safety joint
- Desiccators with flat flange and MOBILEX screw thread joint

Due to precision manufacture of DURAN desiccators of the same type and size spare parts such as bases, lids, stopcocks and closures are fully interchangeable. By combination of different parts the DURAN system offers different variations, according to individual needs.

These variations are also possible if you wish to combine existing SBW desiccator parts with DURAN parts.

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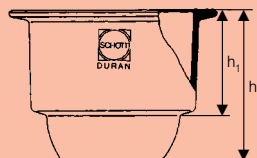
Desiccator bases

with flat flange, for all types of lids,
DIN 12 491

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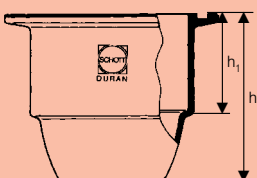
Desiccator bases

with ring-grooved flange,
for all types of lids

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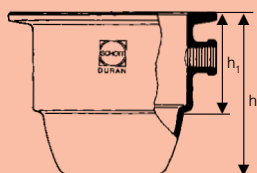
Desiccator bases

with plane flange, with NS
side outlet, NOVUS type

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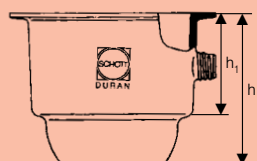
Desiccator bases

with flat flange, screw thread outlet,
MOBILEX type

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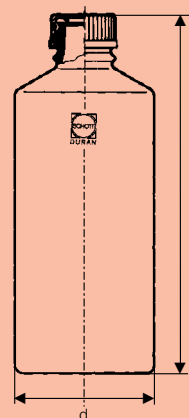
Roller bottles for cell cultures
with DIN-thread
with plastic screw-cap and pouring
ring in polypropylene, blue

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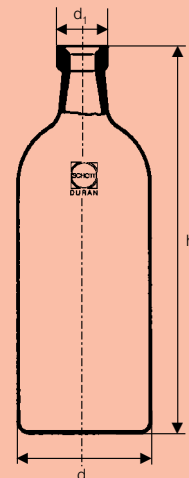
Bottles with rolled flange
without closure

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Bottles with rolled flange
with clamp closure

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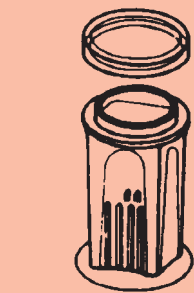
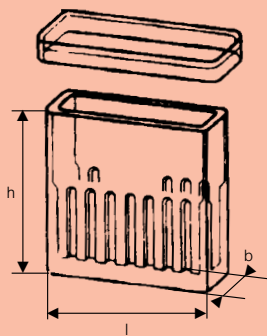
Staining jar
for 10 microscope slides 76 x 26 mm
made of soda-lime glass*

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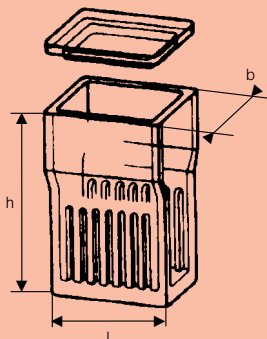
Staining dish, Hellendahl type
for 16 microscope slides 76 x 26 mm
made of soda-lime glass*

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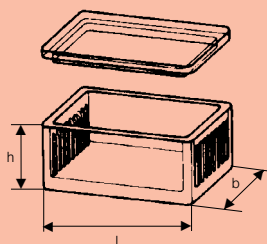
Staining dish, Hellendahl type
for 16 microscope slides 76 x 26 mm
with wide upper section
made of soda-lime glass*

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**Staining dish,
Schiefferdecker type**
for 10 microscope slides 76 x 26 mm
made of soda-lime glass*

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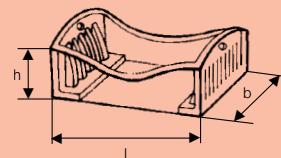
Staining tray, DURAN

for 10 microscope slides 76 x 26 mm
or any width up to 52 mm

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Glass box

with cover for staining tray
made of soda-lime glass*

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DURAN Volumetric Flasks

Volumetric flasks which are used to measure accurately specific liquid volumes, are, like most other volumetric glassware, indispensable aids to quantitative analysis. The chemist uses them principally in the preparation and storage of standard solutions. DURAN volumetric flasks are made of the chemically resistant borosilicate glass 3.3 manufactured by Schott.

They are calibrated to contain ("In") at the reference temperature of 20 °C. The tolerances of volume content of class A volumetric flasks correspond to the accuracy limits of Federal German weights and measures requirements as well as to recommendations of ISO¹ and DIN².

¹ International Organization for Standardization

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Volumetric flasks

with one graduation mark and stopper made of polyethylene
ISO 1042, DIN 12664, class A

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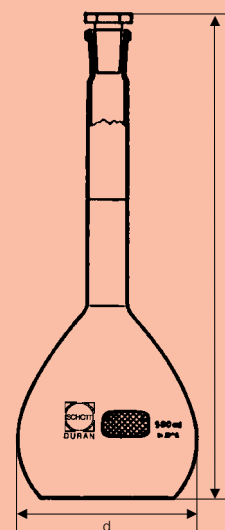
DURAN Measuring and Mixing Cylinders

Measuring cylinders serve simultaneously to receive and measure different amounts of liquid. Mixing cylinders are used when diluting solutions and when mixing a number of components in pre-determined volume ratios. DURAN measuring and mixing cylinders are made of borosilicate glass 3.3 and are therefore very resistant both to mechanical stresses and thermal shock. The large hexagonal base with **embossed under-surface** increases stability and prevents the cylinder

rolling when in the horizontal position. Cylinders have the same wall thickness over the whole of their graduated portion. This eliminates errors due to taper. They are calibrated to contain (“In”) at the reference temperature of 20 °C. Accuracy limits of measuring and mixing cylinders have been fixed according to DIN¹ 12 680, DIN 12 685 and ISO² 4788.

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² International Organization for Standardization

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Mixing cylinders with hexagonal base

with graduation
with standard ground joint
and plastic stopper
DIN 12685, ISO 4788

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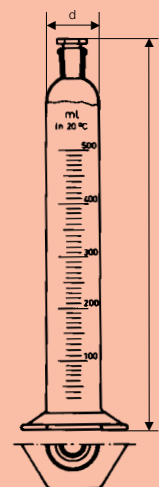
Measuring cylinders with hexagonal base

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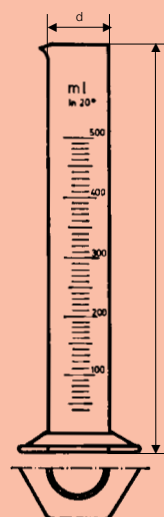
**Zylindermensuren niedrige Form**

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Measuring cylinders, graduated low form

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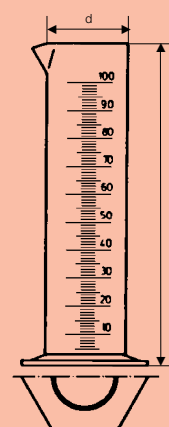
DURAN-Burettes

Burettes are used exclusively for titration. The unknown volume of liquid required in the titration can be read off accurately when the reaction end point has been achieved. DURAN-Burettes are made of the chemically resistant borosilicate glass 3.3 by Schott. They are calibrated to deliver („Ex“) at the reference temperature of 20 °C. The tolerances of burettes on volume content correspond to the accuracy limits of ISO and DIN. The accuracy limits for Schott class B types are about 1½ times as wide as for

class AS. This means they are better than required by ISO or DIN. By specifying a class “AS” within the framework of their 15th modifications order, the Federal German weights and measures authorities have recognized that an overwhelming portion of volumetric measurements, particularly in clinical laboratories, is carried out on water or on dilute aqueous solutions. Thus, apparatus with the same accuracy limits, but with run-out times considerably shorter than hitherto demanded, are now admitted for calibration.

Inhalt Capacity	Fehlergrenze Klasse AS Accuracy limits class AS suitable for official calibration	Fehlergrenze Klasse B Accuracy limits class B	
	DIN 12 700	ISO 385 DIN 12 700	Schott
ml	± ml	± ml	± ml
1	0,01	–	–
2	0,01	–	–
5	0,01	–	–
10	0,02	0,05	0,03
25	0,03	0,05	0,04
50	0,05	0,1	0,08
100 ¹	0,08	0,2	0,15

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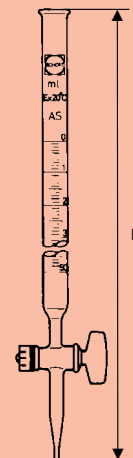
Burettes

with straight stopcock, with
Schellbach stripe, main graduations
encircle the tube, class AS,
DIN 12700
waiting time 30 seconds

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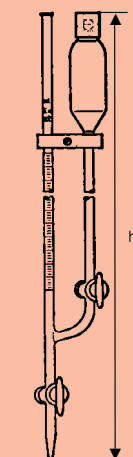
Automatic burettes

according to Pellet, with side stop-
cock, with Schellbach stripe, main
graduations encircle the tube,
class AS, DIN 12 700
waiting time 30 seconds
with reservoir bottle 2,000 ml and
blowball

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DURAN-Bürettenflaschen

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Reservoir bottle (Burette bottle)

with standard ground,
Neck socket size 29/32, DIN 12463

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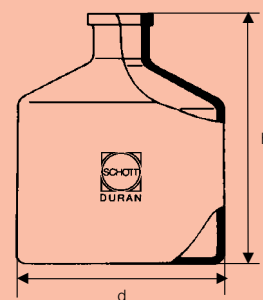
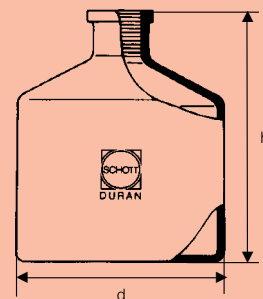
Pipettes

Pipettes are used for accurate measurement and transfer of liquids. Bulb pipettes will only deliver specific quantities.

Graduated pipettes allow take-up of different liquid volumes and delivery of the same or different quantities. Graduated and bulb pipettes are AR-Glas products.

They are calibrated to deliver („Ex“) at the reference temperature of 20 °C. In the case of calibrated pipettes and class AS pipettes, the tolerances on volume content correspond to the accuracy limits of ISO and DIN. The accuracy limits for Schott class B

types are about 1½ times as wide as for class AS. This means that they are better than required by ISO or DIN. By specifying a class „AS“ within the framework of their 15th modifications order, the Federal German weights and measures authorities have recognized that an overwhelming portion of volumetric measurements, particularly in clinical laboratories, is carried out on water or on dilute aqueous solutions. Thus, apparatus with the same accuracy limits, but with run-out time considerably shorter than hitherto demanded, are now admitted for calibration.

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**Tolerances for graduated
measuring pipettes calibrated
to deliver**

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**AR-Glass measuring pipettes
for partial outflow**
classe B
ISO 835, DIN 12 695
with line divisions
graduated with zero at the top

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Klasse B

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AR-Glass measuring pipettes for complete outflow

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Klasse AS, DIN 12 697, ISO 835

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Wartezeit 15 s

AR-Glass measuring pipettes for complete and partial outflow

class AS, DIN 12 697, ISO 835

main graduations encircle the tube

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Screwfilters with interchangeable filter discs

With just 3 filter sizes, each having 4 filter discs of varying porosity, it is possible to obtain 12 different filter rates. DURAN screwfilters have many advantages over conventional filter units:

- interchangeable filter discs
- safe and simple removal of the filtered material
- longer lasting discs, as no damage is caused by scraping off the filtered material

- easy cleaning of the filter disc from both sides
- slit sieve (**21 340 31**) can be used to support membrane- and paper filters
- space saving
- economical; both filter discs and units can be ordered individually, as required

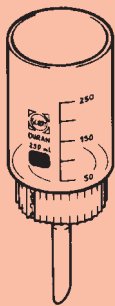
Important:
Place the filter disc between 2 FKM seals.

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Filter head with PP¹ funnel and 2 FKM¹ seals

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² Suitable slit sieves **21 340 31** as support for membrane and paper filters.

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Filter discs with glass rim must be ordered separately.

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Filter disc with glass rim patented

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Fig.: Absorption efficiency of two gas wash bottles
A without gas filter and
B with gas filter plate

DURAN®- gas wash bottles

If the distribution of the gas in the liquid in gas wash bottles is effected by means of a filter plate, the absorption efficiency can be significantly increased. Gas wash bottles of this

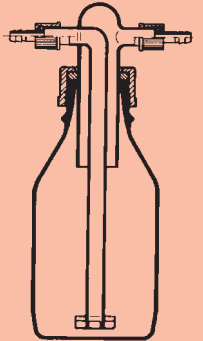
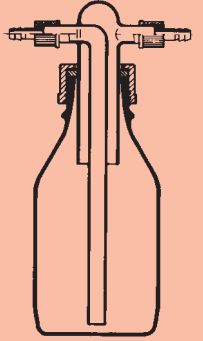
type still operate reliably even at high flow rates. The diagram shows the efficiency of a gas wash bottle with and without glass filter plate.

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Gas washing bottle

screw-cap system
with Drechsel-head
without filter disc
adjustable immersion

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Gas washing bottle

screw-cap system
with Drechsel-head
with filter disc
adjustable immersion

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Filtering flasks and bottles with side-arm socket¹

for vacuum use

DIN 12 476, ISO 6556

These filtration flasks have a ground socket 17,5/26 to accommodate vacuum tubing of 15 to 18 O. D. (e. g. 6 x 5 mm or 8 x 5 mm, DIN 12 865).

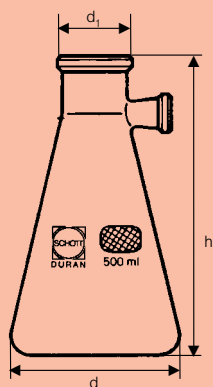
Filtration flasks with socket have not only made work in preparation and analytical laboratories easier and simpler, but have also reduced accident risk.

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Connecting to vacuum

Appropriate vacuum tubing is inserted into the side-arm of the filtration flask until it is firmly held by the socket's conical shape. After connecting the

tubing to a water pump, its inserted end is pulled more tightly into the socket and seals off the connection.

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Filtering flasks, Erlenmeyer shape, with glass hose connection¹, for vacuum use, ISO 6556

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Filtering flasks¹, plastic-coated (DURAN Protect)², with plastic hose connection and plastic tubulature, complete³
for vacuum use

The plastic coating provides protection against mechanical damage and in the event of implosion will hold the broken glass pieces together. The copolymer based plastic material remains transparent and adheres to the glass surface.

The bottles can be cleaned in conventional laboratory apparatus washing machines.

Autoclavable up to 135 °C steam; do not use distilled water or vacuum for cooling.

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¹ These filtering bottles comply with the regulations of the German Gerätesicherheitsgesetz.

² Thermal and chemical durability of VMQ see page 180.

³ With these bottles we supply an assembly set giving you the choice of using either a plastic hose connection or a plastic tubulature.

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DURAN flat flange range by SCHOTT

For many years laboratories of all types have valued DURAN flat flange reaction vessels for their multi-purpose use.

SCHOTT supplies an extensive range of blanks and finished items which can provide the optimum solution for each individual application, whether it be reaction, distillation, evaporation or desiccation.

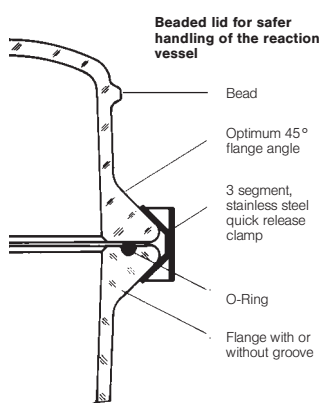
SCHOTT's flat flange and reaction vessels have a number of special features:

- the high angle of taper (45°) results in a very robust design

- the 3 segment, stainless steel quick release clamp ensures easy and safe handling
- the body and lid of the flat flange and reaction vessels have the same nominal diameter, making them interchangeable
- proven flange design – flat ground
 - with or without groove, suitable for any seal required
- beaded lid design for safer handling: vessels can be changed without taking the apparatus apart.

Important:

All assemblies are resistant to vacuum and approved for positive operating pressure as mentioned.

Round-bottom jacketed vessels

SCHOTT supplies round bottom cylinders for the manufacture of jacketed vessels. The matching inner and outer vessels are available in a standard length of 400 mm in the various diameters.

They can be cracked off at the appropriate length so that reaction vessels can be manufactured with capacities ranging from 250 ml to 10 l.

Advantages:

- the use of borosilicate glass type 3.3 ensures compatibility in welding components together
- the vessel blank has been designed to ensure that it can be used under pressure or vacuum (as mentioned)
- the outer and inner vessels are carefully matched to ensure that the flow distribution of the heating fluid in the jacket is uniform.

Notes for the further processor:

Jacketed vessels, flat flange reaction vessels and cover blanks

Tips for apparatus makers

1. The figures given in the operating pressure tables are only valid for Schott's original vessel blanks, on condition that no subsequent reduction in wall thickness occurs.
2. For glass articles which differ from the original shape after further processing, the approved operating data/stress data must be determined in accordance with the technical rules and regulations.

Accessories:

For flat flange reaction vessels, two sealing possibilities are offered:

- a. O rings for operations under superatmospheric pressure and vacuum **up to 200 °C**
 - easy to open

- the cover does not seize, not even after extensive operations under vacuum and higher temperatures
- the supporting surfaces do not have to be greased.

- b. Grease for operations under superatmospheric pressure and vacuum **above 200 °C**
 - the groove acts in this case as an ideal greasing groove
 - no contamination of the substance.

The new stainless steel quick seals with three holding segments provide even distribution of the contact pressure by optimum adaptation.

The chromium-nickel-steel holder comprising two tensioning rods is intended for incorporating the reaction vessels or the covers, for example, in fixing walls. If, for example, the cover or vessel must be exchanged, this can be done without dismantling the entire apparatus.

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Shape-retentive O-rings

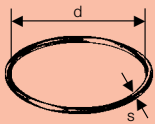
FEP – seamlessly coated elastomer O-rings – with silicone core

These O-rings from SCHOTT consist of a rubber-elastic core and a FEP coating which encloses the rings seamlessly. The combination of these high-quality materials ensures the elasticity of a conventional O-ring. The chemical resistance of FEP (Tetrafluoroethylene-hexafluorpropylene/copolymer FEP) is equal to PTFE, i. e. it is resistant to almost all chemicals.

FEP's temperature resistance ranges from - 50 °C to + 200 °C.

Advantages of the O-rings:

- excellent recovery through elastomer core making O-rings reusable
- excellent chemical resistance, compatible with most liquids/chemicals
- temperature-stable
- solvent-resistant
- physiologically completely harmless

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Dropping funnels, cylindrical, graduated

with pressure equalisation tube, with standard ground cone, with standardized hollow key and key retaining device, with standard ground glass stopper, DIN 12567, ISO 4800

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Separating funnels, conical

with standardized hollow key and key retaining device, with plastic stopper, DIN 12451, ISO 4800

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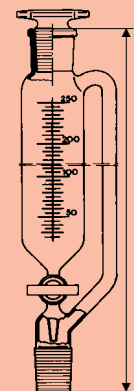
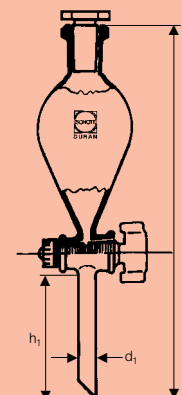
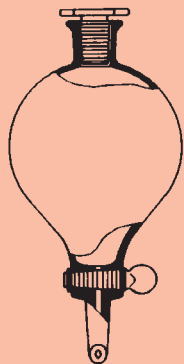
Separating funnels

with NS stopcock and octagonal plate NS stopper, DIN 12450

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KPG stirrer bearings

interchangeable, bearing surface
ground, with standard ground cone

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KPG stirrer bearings

interchangeable,
bearing surface ground

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KPG stirrer bearings

interchangeable, bearing surface
ground, with screwthread and
standard ground cone

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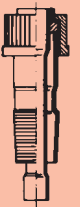
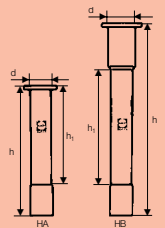
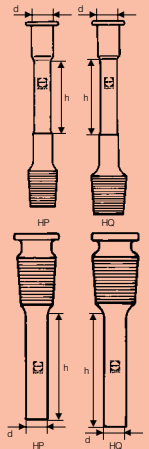
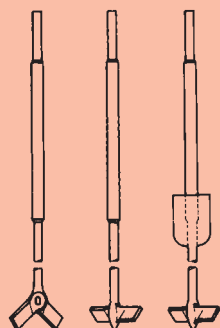
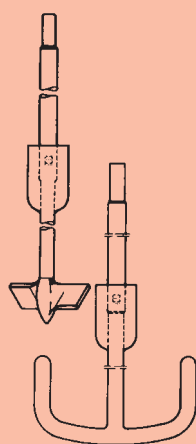
KPG stirrer shafts, ø 10 mm
interchangeable
bearing surface ground and polished

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KPG stirrer shafts, ø 16 mm
interchangeable, bearing surface
ground and polished (particularly
suitable for flat-flange reaction
vessels)

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KPG stirrer shafts
interchangeable
bearing surface ground and
polished

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Distillation thermometers

(solid stem thermometers)
calibrated totally immersed

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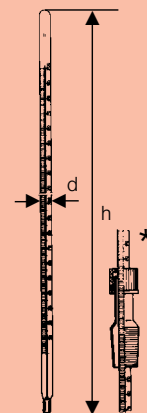
* Additional required screw-caps, seals and screw thread tubes with standard ground cone
see page 113-115.

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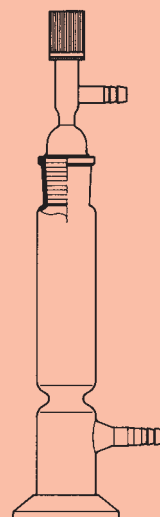
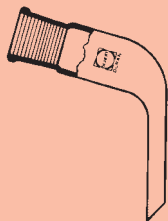
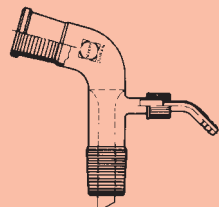
Receiver adapters, plain, bent with standard ground socket

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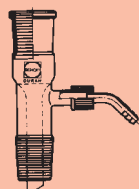
Receiver adapters, bent with vacuum connection, with 2 standard ground joints, with screw- on plastic hose connection

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Receiver adapters, straight with vacuum connection, with 2 standard ground joints, with screw- on plastic hose connection

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Normschliff-Hülsen

ISO 383, DIN 12249

Standardized conical joint-sockets

ISO 383, DIN 12249

Joints femelles à rodage standard

ISO 383, DIN 12249

Hembras esmeriladas

ISO 383, DIN 12249

Best.-Nr. Cat. No.	Nenngröße NS Size designation	Schliffkegel d ₁ h mm mm	Rohr Tube d	l mm	Originalpckg./Stück Stand. pack/quantity	
24 111 01	5/13	5	13	8	115	10
24 111 02	7/16	7,5	16	10	115	10
24 111 03	10/19	10	19	13	115	10
24 111 04	12/21	12,5	21	15	115	10
24 111 06	14/23	14,5	23	17	115	10
24 111 07	19/26	18,8	26	22	125	10
24 111 08	24/29	24	29	28	135	10
24 111 09	29/32	29,2	32	33	135	10
24 111 11	34/35	34,5	35	40	150	1
24 111 12	45/40	45	40	50	150	1

Normschliff-Kerne

ISO 383, DIN 12249

Standardized conical joint-cones

ISO 383, DIN 12249

Joints mâles à rodage standard

ISO 383, DIN 12249

Machos esmerilados

ISO 383, DIN 12249

Best.-Nr. Cat. No.	Nenngröße NS Size designation	Schliffkegel d ₁ h mm mm	Rohr Tube d	l mm	Originalpckg./Stück Stand. pack/quantity	
24 113 01	5/13	5	13	4	120	10
24 113 02	7/16	7,5	16	6	120	10
24 113 03	10/19	10	19	8	120	10
24 113 04	12/21	12,5	21	11	120	10
24 113 06	14/23	14,5	23	13	120	10
24 113 07	19/26	18,8	26	17	150	10
24 113 08	24/29	24	29	22	150	10
24 113 09	29/32	29,2	32	26	150	10
24 113 11	34/35	34,5	35	30	150	1
24 113 12	45/40	45	40	40	150	1

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DIN 12257

Adapters, reduction and expansion

with standard ground conical joints
DIN 12257

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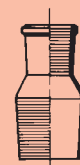
Clips for conical joints of nichrome wire

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Keck assortment for clips for conical joints, POM¹

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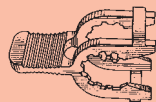
Keck assortment for clips for conical joints, metal

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Keck assortment for clips for spherical joints, POM¹

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Glass flat-head stoppers

with standard ground cone,
octagonal, DIN 12252

massiv/solid/massifs/macizo			
21 624 03	–	12/21	10
21 624 04	–	12/21	10
21 624 06	24 624 06	14/23	10
21 624 07	24 624 07	19/26	10
–	24 624 08	24/29	10

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Glass flat-head stoppers

with short standard ground joint,
octagonal, DIN 12252

massiv/solid/massifs/macizo			
21 625 06	24 625 06	14/15	10
21 625 08	24 625 08	24/20	10

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DURAN glass stoppers*

ground conical, for reagent bottles
(oxygen bottles according to
Winkler)

21 622 06	24 622 06	50	14/23	10
21 622 07	24 622 07	100	19/26	10

* Passend für Standflaschen s. S. 28-31.

* Suitable for reagent bottles see page 28-31.

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Important note: When DURAN glass stoppers are used with soda-lime glass reagent bottles, they may become stuck in the neck if temperature differences exceed 30 °C.

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Deckelstopfen, braun

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Glass flat-head stoppers, amber

with standard ground joint,
octagonal, DIN 12252

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Plastic screw-caps with aperture, red

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SVS screwthread adapter couplings for 2 x GL

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Best.-Nr. Cat. No.	Nenngröße für Nominal size for	Originalpckg./Stück Stand. pack/quantity
29 226 055	2 x GL 14	1
29 226 065	2 x GL 18	1
29 226 095	2 x GL 25	1
29 226 085	2 x GL 32	1

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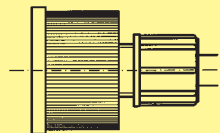
Sealing rings (silicone + PTFE) for polished spherical joints

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Silicone rubber sealing for GL threads with PTFE washer

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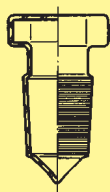
Screwthread tubes with cone, GL/cone

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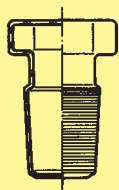
Hexagonal stoppers, pointed, DG ground

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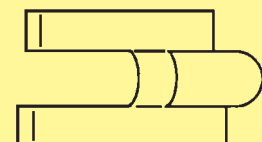
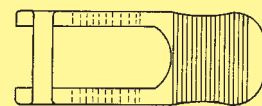
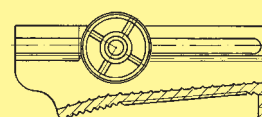
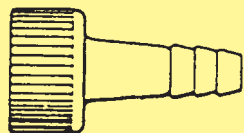
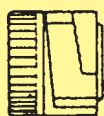
KECK tubing part for adapters KA, with EPDM sealing 16 mm

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KECK screw cap for adapter KA, with sealing EPDM 12 mm

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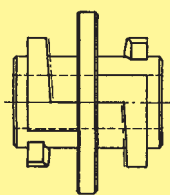
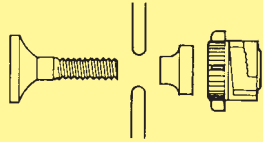
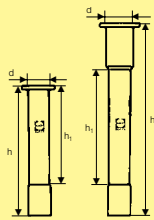
**CPG tubes, ends cut,
length 500 mm, shortest length
to use 250 mm**

**Tubes KPG, extrémités coupe
scie diamantée, long. 500 mm,
long. utile 250 mm**

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CPG stirrer bearings, interchangeable

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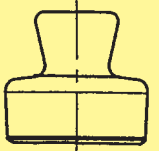
Stopcock side arms, with extension jet

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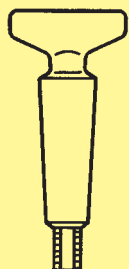
Stopcock keys threaded, SBW glass, without bore, unground

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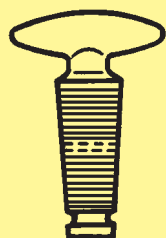
Stopcock keys threaded, SBW glass, long for double bore, without bore, unground

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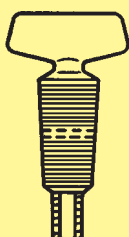
Stopcock keys threaded, SBW glass, ground

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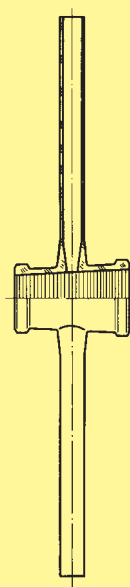
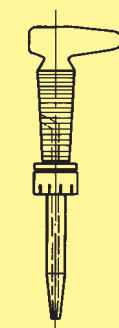
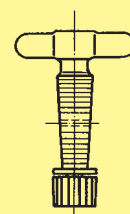
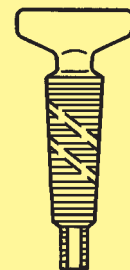
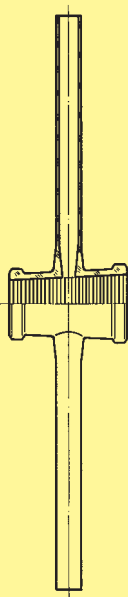
Single way stopcocks, ground, without keys

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Single way stopcocks, ground and polished, without keys

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Single way stopcocks, complete with DURAN key with groove

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Single way stopcocks, complete with SBW key

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Single way stopcocks, taper 1 : 5, with PTFE keys

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Stopcocks, two way, double oblique bore, ground, without keys

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Stopcocks, two way, double oblique bore, ground and polished, without keys

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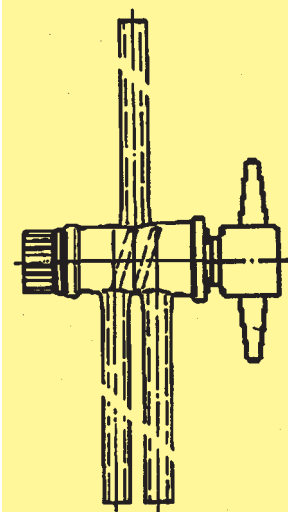
Stopcocks, two way, double oblique bore, complete with SBW-keys

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**Stopcocks, two way,
double oblique bore,
complete with PTFE-keys**

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**3-way stopcocks, complete
with SBW-keys, T-bore**

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**3-way stopcocks, complete
with PTFE-keys, T-bore**

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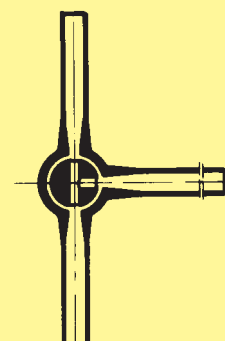
**Capillary 3-way stopcocks,
DUROBAX, complete with
SBW-keys, T-bore**

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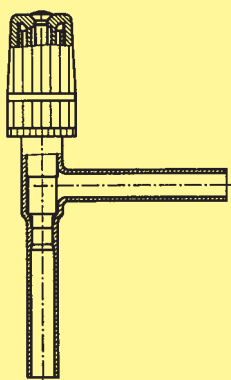
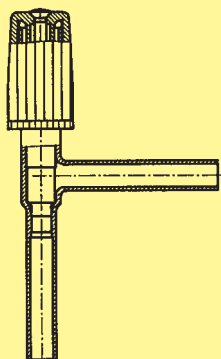
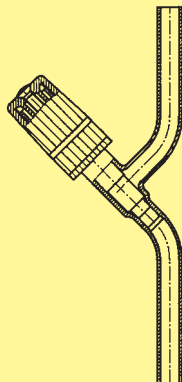
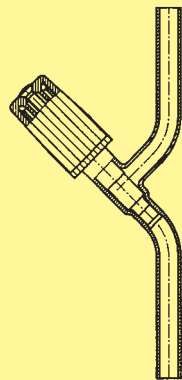
**Burette stopcocks, straight,
AR glass, ground, without keys**

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**Burette stopcocks, straight,
AR glass, ground and polished,
without keys**

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**Burette stopcocks, straight,
fine ground, without keys**

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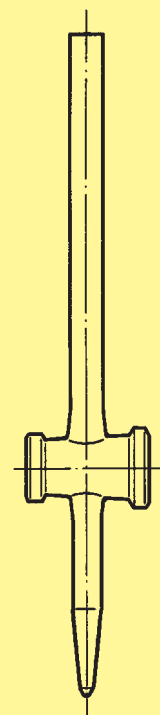
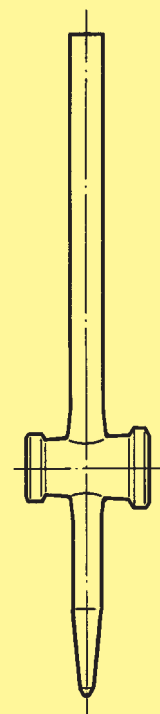
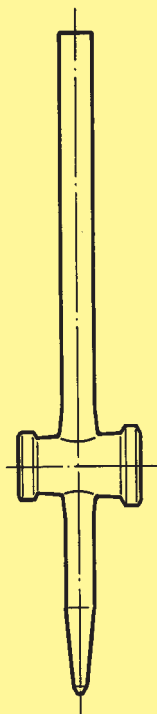
**Burette stopcocks, straight,
ground, without keys**

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**Burette stopcocks, straight,
ground and polished,
without keys**

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**Burette stopcocks, straight,
AR glass,
complete with SBW-keys**

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**Burette stopcocks, straight,
AR glass,
complete with PTFE-keys**

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**Burette stopcocks, straight,
complete with SBW-keys**

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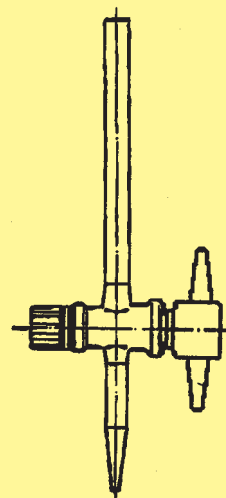
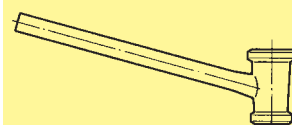
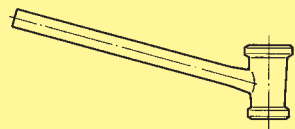
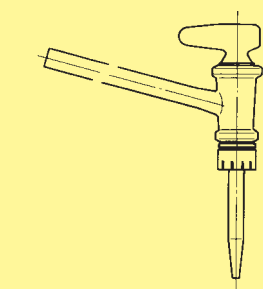
**Burette stopcocks, angled,
ground and polished,
without keys**

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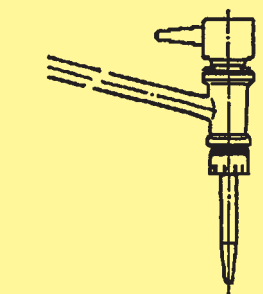
**Burette stopcocks, angled,
AR glass,
complete with SBW-keys**

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**Burette stopcocks, angled,
AR glass,
complete with PTFE-keys**

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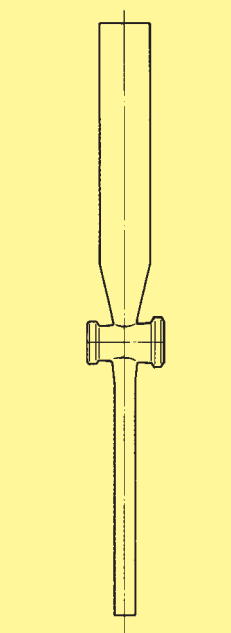
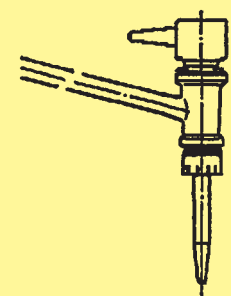
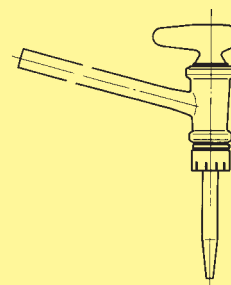
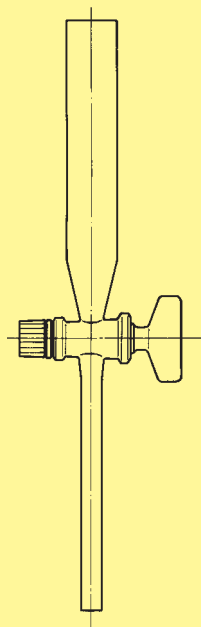
Separating funnel stopcocks, Gilson, with SBW-keys

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Separating funnel stopcocks, Gilson, with PTFE-keys

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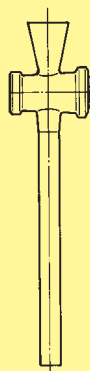
Separating funnel stopcocks, with short inlet, ground, without keys

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Separating funnel stopcocks, with short inlet, ground and polished, without keys

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Separating funnel stopcocks, with short inlet, with SBW-keys

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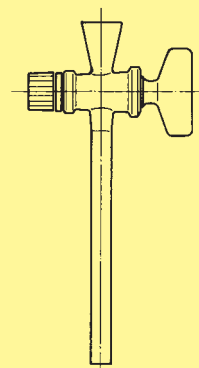
Separating funnel stopcocks, with short inlet, with PTFE-keys

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Gas washing bottle lower part, Drechsel type

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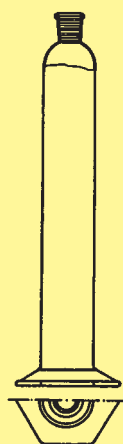
with standard ground joint, without stopper, without graduation, for DIN 12685, ISO 4788

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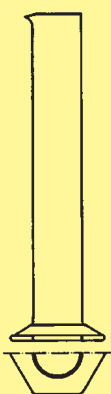
Measuring cylinder

without base, tall form, with spout,
flat bottom, without graduation

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Measuring cylinder

low form

with hexagonal base, with spout,
without graduation

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**DURAN
Laboratory glassware
Edition I/1997**

**All items are subject to technical
and standards related changes.**

In cases of changes in DIN specifications, due, for example, to assimilation to ISO recommendations, our dimensions will be correspondingly adjusted over a reasonable time period.

Delivery through authorized dealers.

We feel unable to grant any rebates on orders of a net value below DM 500,-.

Articles for which standard pack quantities are given, can be supplied only in the quantities mentioned,

these quantities being at the same time the minimum number of pieces to be purchased. Exceptions to the above are specially marked.

Please make your arrangements accordingly.

The terms of delivery and payment are always applicable in their latest edition.

Reproduction of this catalogue, in whole or in part, is permitted only with our written consent.

As from the date this catalogue takes effect, all issues published previously will be void.

Trademarks

SCHOTT, CERAN, DURAN, DUROPLAN, FIOLAX, KPG, SUPREMAX, PRODURAN, SBW, STERIPLAN and WERTEX are registered trademarks of SCHOTT GLASWERKE, Mainz.

AR-GLASS (soda-lime) is a registered trademark of SCHOTT-ROHRGLAS GmbH, Bayreuth.

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DURAN by Schott

DURAN is a borosilicate glass type 3.3 as specified in international standard ISO 3585 and DURAN products meet the most important international standards, for example the German¹, British, American and French standards. Maximum chemical resistance, minimum thermal expansion and the resulting high resistance to thermal shock are its principal properties. This optimum physical and chemical

performance makes DURAN the ideal material for use in the laboratory and for large-scale industrial plants in the manufacture of chemical plant. Moreover, it is widely used on an industrial scale in all other areas of application in which extreme heat resistance, resistance to thermal shock, mechanical strength and exceptional chemical resistance are required.

Definition

Glass is an inorganic mixture fused at high temperature which solidifies on cooling but does not crystallize. Its basic components, network formers and modifiers are present in the form of oxides in ordinary glasses. Typical glass formers (network formers) are silica (SiO_2), boric acid (B_2O_3), phosphoric acid (P_2O_5) and, under certain circumstances, also aluminium oxide (Al_2O_3).

These substances are capable of absorbing (dissolving) certain amounts of metal oxides without losing their glassy character. This means that the incorporated oxides do not participate as glass formers but modify certain physical properties of the glass structure as „network modifiers“.

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¹ DIN and EN standards can be purchased from: Beuth Verlag GmbH, Postfach 11 45, D-10772 Berlin
ISO denomination according to ISO 4791/1

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Glass structure

A large number of chemical substances solidify from the liquid state of the melt in the form of a glass. The glass formation depends on the cooling rate and requires mixed types of bonding (covalent bond and ionic bond) between the atoms or atom groups. As a result, glass-forming products have a strong tendency while still in the melt for three-dimensional crosslinking by polymerisation in a largely random manner. Crystals are formed when the individual atoms form a regular three-dimensional arrangement in what is known as a

“crystal lattice”, as soon as the particular substance changes from the liquid to the solid state. Glass, however, on cooling from the liquid state, forms a largely spatially random “network”. The main components which participate in the glass formation are therefore called “network formers”. Ions can be incorporated in this network of glass-forming molecules, as a result of which they tear up the network in certain places and modify the network structure and thus glass properties. That is why they are called “network modifiers”.

Chemical composition of DURAN®

DURAN glass which is used in the laboratory and in industrial scale chemical plant, due to its excellent chemical and physical properties, has the following typical composition:

81% by weight SiO_2
13% by weight B_2O_3
4% by weight $\text{Na}_2\text{O}/\text{K}_2\text{O}$
2% by weight Al_2O_3

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Chemical properties

The chemical resistance of glass is more extensive than that of all other known materials. Borosilicate glass is highly resistant to water, acids, salt solutions, organic substances and also halogens such as chlorine and bromine. It also has relatively good resistance to alkaline solutions. Only hydrofluoric acid, concentrated phosphoric acid and strong alkaline solutions cause appreciable surface removal of the glass at higher temperatures.

Hydrolytic resistance according to ISO 719 (98 °C)

Hydrolytic resistance by the glass grain method class ISO 719-HGB 1 (corresponds to the previous DIN 12111, hydrolytic class 1).

Hydrolytic resistance according to ISO 720 (121 °C)

Hydrolytic resistance by the glass grain method class ISO 720-HGA 1.

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Fig. 2:
Hydrolytic attack on DURAN as a
function of time (100 °C) approx. to
ISO 719.

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**Acid resistance according to
DIN 12 116**

Acid class 1

**Acid resistance according to
DIN-ISO 1776**

Removal of $\text{Na}_2\text{O} < 100 \mu\text{g}/\text{dm}^2$
The attack curves in Figure 3 show a maximum in acid ranges between 4-7 N (HCl, for example in the azeotrope containing 20.2% by weight). At higher concentrations, the reaction rate decreases significantly, so that the layer thicknesses which are

attacked are only in the range of a few thousandths of a millimetre after years. Thus, the mechanisms of acid attack are not relevant given the wall thicknesses used in practice. Therefore, it is rightly called “acid-resistant borosilicate glass”.

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Fig. 3:
Acid attack on borosilicate glass
as a function of acid concentration.

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Alkali resistance according to DIN 52 322

Alkali resistance class DIN 52 322-A2 (corresponds to alkali resistance class ISO 695-A2). The surface removal curve in Figure 4 shows that visible attack on the glass surface does not take place until above about 60 °C. At lower temperatures (incorporation of coolers in the NaOH cycle), the reaction rates are so low that hardly any reduction of the wall

thickness takes place over a period of years. Long-term tests have shown that the use of NaOH having a concentration of 1 mol/l (corresponds to 4% strength by weight sodium hydroxide solution, pH 14) at an operating temperature of 50 °C gives a glass surface removal of 1 mm after 25 years in a continuous flow through a glass pipeline.

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Fig. 4:
Alkali attack on DURAN as a function of temperature.

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Ecologically harmless

Special glass is considered completely safe ecologically. Glass is manufactured from natural raw materials (sand, calcium carbonate and sodium carbonate) and does not

contain any materials which could be released and constitute a hazard to man or the environment. Glass can be reused several times and disposed of without difficulties.

Use in microwave apparatus

DURAN laboratory glass is suitable for use in microwave apparatus.

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Sealing DURAN to related glasses

With its heat expansion of $3.3 \times 10^{-6} \text{ K}^{-1}$, DURAN is unsuitable for direct seals with sealing metals. For seals with tungsten or tungsten glasses, the intermediate sealing glass 8448 must

be used. For seals with glasses of the molybdenum/Kovar range, two intermediate sealing glasses as shown in the figure below are recommended.

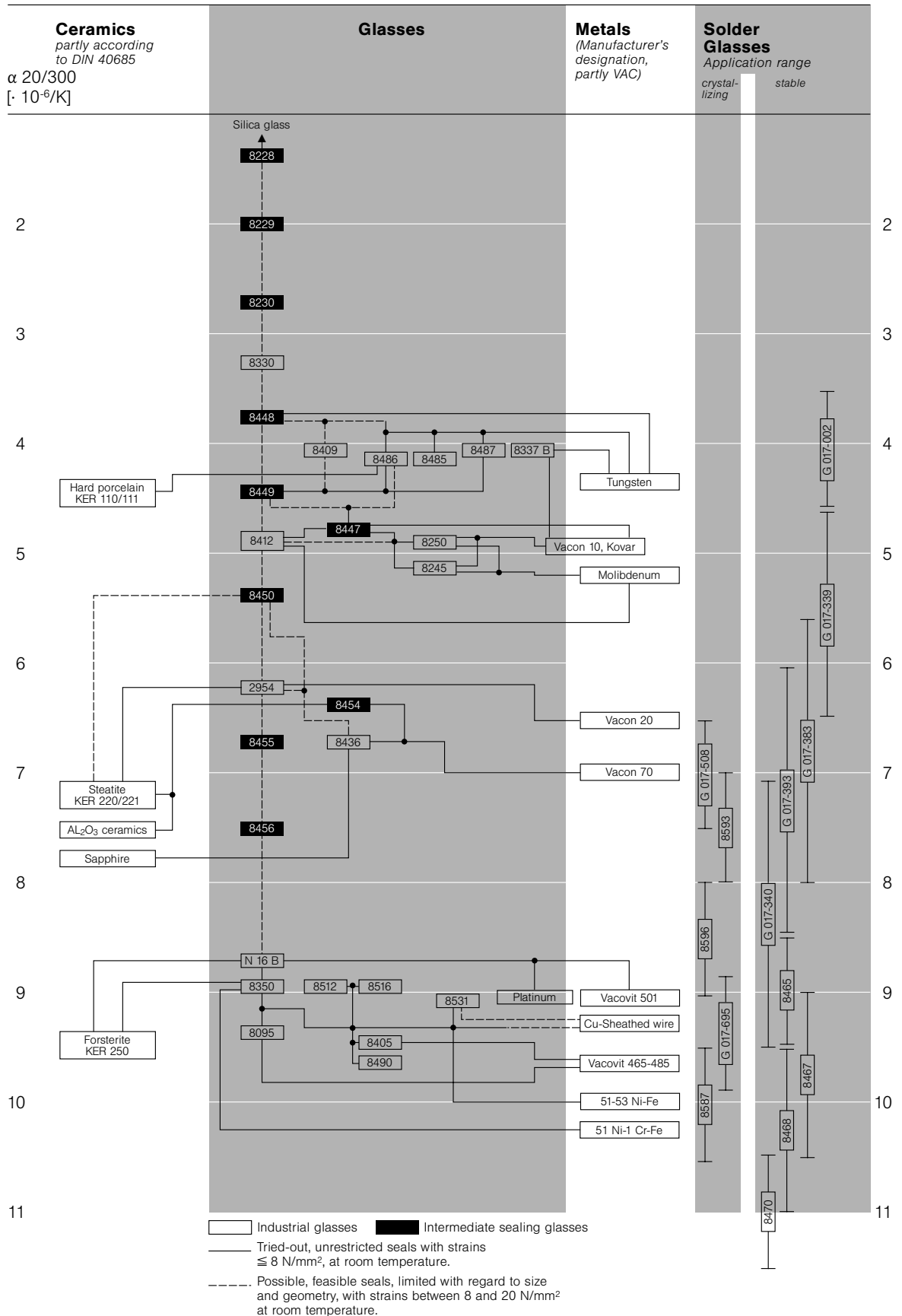
Fig. 5: Recommended technical glass seal combinations with other glasses, important metals, alloys and ceramics.

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Optical properties

Spectral range in which the absorption of DURAN is negligibly small:
approx. 310 to 2,200 nm absorption of brown glass up to about 500 nm

SCHOTT laboratory glasses do not show any significant absorption in the visible spectral range. Thus, DURAN appears transparent and colourless. Fairly large layer thicknesses (axial view through pipes) appear greenish. When working with light-sensitive substances, the surfaces can be coloured brown.

The result is strong absorption in the short-wave region. In photochemical processes, light transmission of DURAN in the ultraviolet region is of particular importance. The degree of transmission in the UV region shows that photochemical reactions can be carried out, for example chlorinations and sulphochlorinations. The chlorine molecule absorbs in the range from 280 to 400 nm and thus serves as a carrier of the energy of radiation.

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Physical properties and characteristics for the processing of DURAN by glassblowers

Maximum use temperature: 500 °C
Transformation temperature DIN-ISO 3585: 525 °C

Upon cooling from the melting temperature (10¹ to 10³ dPa s), the glasses go through a wide range of viscosity, which ranges from a relatively viscous region (at about 10⁴ to 10⁸ dPa s) to the plastic range (10⁸ to 10¹² dPa s) and up into the known elastic-brittle state (> 10¹³ dPa s). Thus, glass shows a constant increase in viscosity upon cooling and has no solidification point as is the case for crystalline substances. For

the processing of glass, certain sections of the viscosity range are of particular importance. In the transformation range, the elastic-brittle behaviour of the glass changes to a noticeably viscous behaviour with increasing temperature, which significantly changes all physical and chemical properties as a function of temperature. Thus, the temperature region of the transformation range determines the elimination of stress during heating and the onset of stresses during cooling of the glass. The position of the transformation range is indicated by the transformation temperature T_g DIN 52324.

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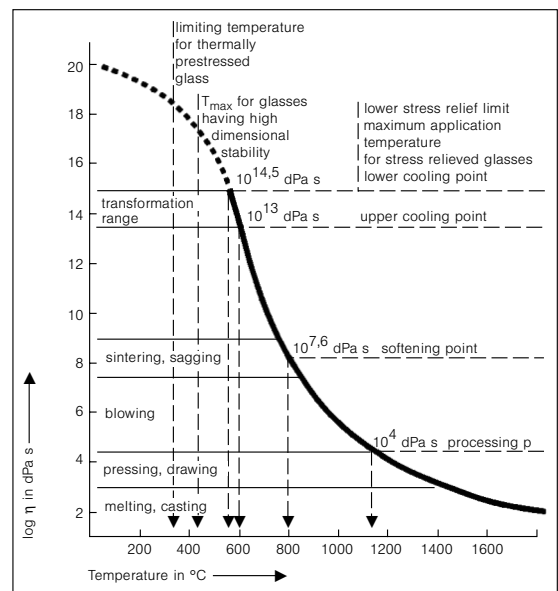
Fig. 7:
 Normal temperature dependence/viscosity curve of, for example, DURAN; viscosity ranges of important processing techniques, position of fixed points of viscosity and various limiting temperatures.

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To eliminate permanent stresses, DURAN is heated to a maximum temperature of 550 °C and kept at this temperature for 30 minutes.

For the cooling which follows, the cooling rates listed in the table below are recommended.

Temperature range	550 to 480 °C	480 to 400 °C	400 to 20 °C
Glass thickness	Cooling rate		
3 mm	12 deg/min	24 deg/min	to 480 deg/min
6 mm	3 deg/min	6 deg/min	to 120 deg/min
12 mm	0.8 deg/min	1.6 deg/min	to 32 deg/min

Table 1

These data are those of today's knowledge and are intended to provide information on our products and their possible applications. Thus, their meaning is not to ensure certain

properties of the products or their suitability for a particular application purpose. We guarantee good quality in accordance with our general conditions of sale.

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Unless stated otherwise, our products are made of DURAN borosilicate glass. Articles made of other glasses are specifically indicated:

Important:

DURAN is a neutral glass of high hydrolytic resistance and belongs to glass type I, according to DAB 10 and the European Pharmacopoeia.

Name	Melt No.	Linear coefficient of expansion α 20/300 °C 10^{-6} K^{-1}	Transformation temperature °C	Density g/cm ³
DURAN BOROFLOAT	8330	3.3	525	2.23
SUPREMAX	8409	4.1	745	2.57
FIOLAX	8412	4.9	565	2.34
AR-GLAS/ soda-lime- silicate glass	8350	9.1	525	2.50
SBW	8326	6.5	555	2.45

Name	Class of chemical durability		
	Water resistance DIN-ISO 719	Acid resistance DIN 12 116	Base resistance DIN 52322/ISO 695
DURAN BOROFLOAT	1	1	2
SUPREMAX	1	4	3
FIOLAX	1	1	2
AR-GLAS/ Soda-lime- silicate glass	3	1	2
SBW	1	1	1

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Fig. 8: Alkali attack on borosilicate glass related to temperature and calculated from the weight losses. NaOH = 1 mol/l. Exposure time: 1 h.

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Fig. 9: Alkali attack on DURAN related to pH value at 100 °C.

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Cleaning

Laboratory glassware can be cleaned manually in a dipping bath or by machine in a laboratory dishwashing machine. A wide range of cleaners and disinfectant cleaners is available for both methods in specialist stores. As contamination during delivery of our laboratory glassware cannot be totally ruled out, may we suggest that the glass is always washed prior to use. In order to take good care of the laboratory apparatuses, they should be cleaned immediately after use at low temperature, a short residence time and low alkalinity. Laboratory apparatuses which have come into contact with infectious substances are first cleaned and then sterilised with hot air or steam. This is the only way to prevent impurities from sticking onto the glass and prevent damage of the glasses caused by any adhering chemicals.

Manual cleaning

Wiping and scouring with a cloth or sponge, each of which are impregnated with a cleaning solution, are

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generally known. Laboratory glasses must never be treated with abrasive scouring agents or sponges, since this would damage the surface. In the dipping bath procedure, the laboratory glasses are, as a rule, placed into the cleaning solution at room temperature for 20 to 30 minutes, then rinsed with tap water and then with distilled water. Do not lengthen the time of exposure or increase the temperature except in the case of persistent impurities. In the case of laboratory glasses, lengthy times of exposure at more than 70 °C should be avoided in alkaline media, since this may result in destruction of the print.

Machine cleaning

The machine cleaning of laboratory glasses in the laboratory dishwasher is a milder treatment than cleaning in a dipping bath. The glasses only come into contact with the cleaning solution during the relatively short washing cycles, when this solution is sprayed onto the glass surface via spray or injector nozzles.

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Important notes for the user!

When working with glass, the limits of this material upon change in temperature and mechanical stress must be heeded and strict safety precautions must be observed:

- Never subject glassware to abrupt changes in temperature. Thus, do not remove it from the drying oven while hot and place it on a cold or, which is worse, a wet laboratory bench. This is true in particular of thickwalled glassware, such as suction flasks or desiccators.
- Provide apparatuses with a support via a suitable stand and set up free of stress. Use, for example, PTFE corrugated bellows to compensate for stresses or vibrations.
- Glass apparatus (e. g. filtering flasks, desiccators, etc.) must be handled with care when under pressure or vacuum.
- To avoid setting up stresses in the glass, glass vessels under vacuum or pressure should not be heated from one side or with an open flame. The maximum permissible pressures quoted in the catalogue should not be exceeded.

Note on complaints

If in spite of our quality control measures you should have a product complaint in relation to DURAN laboratory glass, then please contact your laboratory dealer or us directly.

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- Before evacuating or pressurising glass vessels always check visually that they are not damaged in any way (heavy scratching or other mechanical damage). Damaged items should not be used under vacuum or pressure.
- Never subject glassware to abrupt changes in pressure, for example never aerate evacuated glassware abruptly. Flat-bottomed glassware (for example Erlenmeyer or flat-bottomed flasks) must not be subjected to pressure or vacuum.

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SCHOTT Glaswerke

Kundenservice Laborglas
Hattenbergstrasse 10
D-55122 Mainz
Germany
Telephone +49-6131-664907
Telefax +49-6131-664051
We will deal with your complaint immediately and contact you as quickly as possible.

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Plastics used with laboratory glass

Temperature resistance
up to °C

PE	Polyethylene	- 40 to + 80
PP ²	Polypropylene	- 40 ⁴ to + 140 ⁴
PBT ²	Polybutyleneterephthalat	- 45 to + 200 ³
PTFE ²	Polytetrafluoroethylene	- 200 to + 260
FEP	Tetra-Fluor-Ethylene/Hexafluor-Propylene	- 200 to + 200
ETFE ²	Partially crystalline ethylene/ tetrafluoroethylene copolymer	- 100 to + 180
VMQ ²	Silicone rubber	- 50 to + 230
EPDM	Ethylene/propylene-diene-rubber	- 45 to + 150
PU ¹	Polyurethane	- 20 to + 135
FKM	Fluorinated rubber	- 20 to + 200
POM	Polyoxymethylene	- 40 to + 90

¹ All laboratory glasses coated with PU may only be cleaned while moist in order to avoid any electrostatic charge which may form.

² This plastic types are recommended by the BGA (Bundesgesundheitsamt) (Federal Department of Health) in accordance with the specifications applicable in each case.

³ Changes in colour may occur at temperature stresses above 180 °C.

⁴ For a short time without mechanical stress.

Chemical resistance of plastics

Classes of
substances
+ 20 °C

	PE	PP	PBT	PTFE/FEP	ETFE	VMQ	EPDM	PU	FKM	POM
Alcohols, aliphatic	++	++	++	++	++	+	+	++	-	+
Aldehydes	+	+	++	++	++	+		++		+
Alkaline solutions	++	++	+/-	++	++	-	++	++	-	+
Esters	+	+	+	++	++	-	++	+	-	-
Ethers	-	-	+	++	++	-	-	+	-	+
Hydrocarbons, aliphatic	-	++	++/+	++	++	-	++	++	++	+
Hydrocarbons, aromatic	-	+	++/+	++	++	-	+	++	++	+
Hydrocarbons, halogenated	-	+		++	++	-	+	-	++	+
Ketones	+	+	+/-	++	+	-	++	+	-	+
Acids, dilute or weak	++	++	++	++	++	-	++	++	++	+
Acids, trated or strong	++	++	+	++	++	-	++	+	++	-
Acids, oxidising (oxidising agents)	-	+	-	++	+	-	-	+	+	-

++ = very good resistance
+ = good to limited resistance
- = low resistance

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SCHOTT GLASWERKE, MAINZ

Terms of Delivery and Payment

The following terms of delivery and payment apply to all goods and services supplied by us. Should our customers' terms of purchase differ from these terms, they will only be accepted if confirmed by us in writing. Acceptance of goods supplied constitutes acceptance of these terms.

1. Prices are in German Deutschmarks at the ex works price in effect on the day of delivery. Unless agreed otherwise, invoices are to be paid immediately without deduction of any discount.

If payment is not made by the due date, legal default proceedings may be taken without further notice. We also reserve the right to charge interest at the rate charged by our own bank at the time and the total balance owing will be due for immediate payment, regardless of any other previous arrangement.

For custom-made items we reserve the right to add a reasonable price supplement and to vary the quantity supplied within reasonable limits. Any quantities in excess of the quantity ordered must be accepted. In orders for blanks and glass-to-metal seals, the amount delivered may deviate from the number actually ordered by around 10%.

2. Every effort is made to maintain quoted delivery dates. However, due to the risks and nature of working with glass, these dates are not in any way binding, unless agreed otherwise. Our contractual obligation to our customers is subject to correct and prompt deliveries from our suppliers.

3. The point of delivery is the works where the goods are manufactured. The place for the fulfilment of payments is Mainz. The goods will be despatched through forwarding agents chosen by us. Once the forwarding agents are in receipt of the goods, the risk passes to the buyer.

4. Packing material is not returnable unless otherwise arranged.

5. Every care is taken in the execution of orders. If, in spite of this, the buyer should have any reason for complaint, notification of this must

be within 14 days of the receipt of the goods. Otherwise the goods will be deemed to have been accepted. Warranty claims for industrial equipment or parts thereof must be made within 6 months of supply in the case of those used on a single shift basis and 3 months for multi-shift use. Parts subject to wear and tear are not covered by warranty. No goods should be returned to us without our prior approval.

If, in spite of our care and attention, the goods supplied prove to be faulty, we reserve the right to either repair or replace them. In the case of damages, including consequential damages, the maximum for which we will be liable is the amount on the invoice and, in the case of glass-to-metal seals, only if more than 3% of the batch are rejected as faulty.

6. All drawings, illustrations, dimensions and weights are approximate, unless otherwise specified. The buyer is responsible for ensuring that any working drawings he submits do not infringe the patent rights of any third party: we will not be held responsible for any resulting compensation claims.

7. Any information that we supply may not be reproduced, made available to third parties, or used for any other than the agreed purpose.

8. The buyer shall make such arrangements for storing and identification of the goods as shall be satisfactory to the seller.

Property in the goods shall not pass to the buyer until payment in full has been received by the seller for all goods sold under the contract and for all goods supplied by the seller under any other contract and until that time the buyer shall hold any goods in a fiduciary capacity on the seller's behalf.

The buyer shall not be prevented by the seller's reservation of title from selling or dealing in the ordinary course of business with any of the goods by all such sales or dealings shall be made by the buyer as agent for the seller and the buyer shall not attempt or purport to pledge or transfer the goods in any way as security to any third party until title to them has passed to the buyer in accordance with these conditions.

Any proceeds of sale of goods sold by the buyer pursuant to above conditions shall be kept by the buyer in a separate account.

The seller shall be entitled, at any time after a default has been made by the buyer on any payment by the due date, to require the buyer by written notice to place at the seller's disposal any goods to which title has been reserved by the seller under these conditions any to pay over to the seller any proceeds of sale in respect of goods which have been sold by the agent for the seller. Any action so taken by the seller shall not prejudice the rights of the seller either with respect to the goods concerned or any other goods or with respect to its action for the price.

9. In the case of sales in countries where the right of ownership described in 8 above does not have the same effect as in German law, the goods remain our property until all contractual obligations of the sale have been fulfilled, and all payments have been made. If the right of ownership does not have the same effect as in German law but the reservation of other rights over the goods is permissible, then we have the authority to exercise these rights. The buyer must be prepared to co-operate with any measures we take for the protection of our right of ownership, or the equivalent right to the goods.

10. All legal relationships formed as a result of the sales agreement are in accordance with German law. Mainz is the place of jurisdiction for all parties and all matters, including those concerning bills of exchange. If Schott appears as a plaintiff, it will also be entitled to start legal proceedings at the premises of the buyer.

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